

LAQUA

Quick Manual

COND meter DS-71




HORIBA
Scientific


CODE : GZ0000272696
October, 2011 © 2011 HORIBA, Ltd.

Preparation

① Connect the electrode to the COND meter

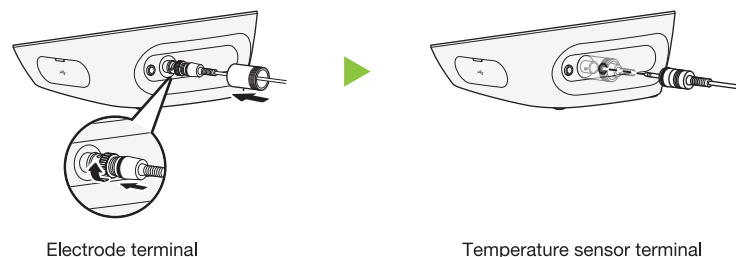


② Press the power button of the COND meter

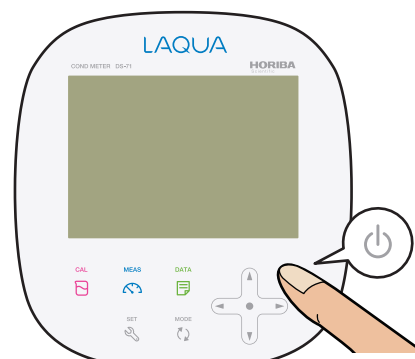

- ☆ Press the power button for 2 seconds.
- ※ Wait for 10 seconds after connecting the AC adapter before pressing the power button.


Set the cell constant

① Enter the set up mode

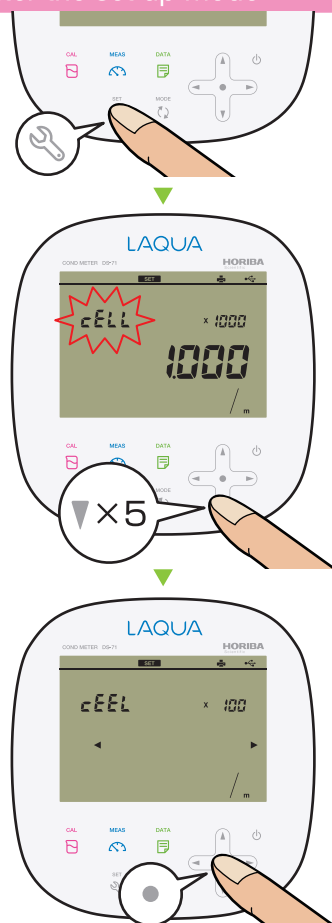

② Set the cell constant


Check the label on the electrode.
Example.) $3552-10D$
 $1.065 \times 100 \text{ m}^{-1} (1\text{cm}^{-1})$

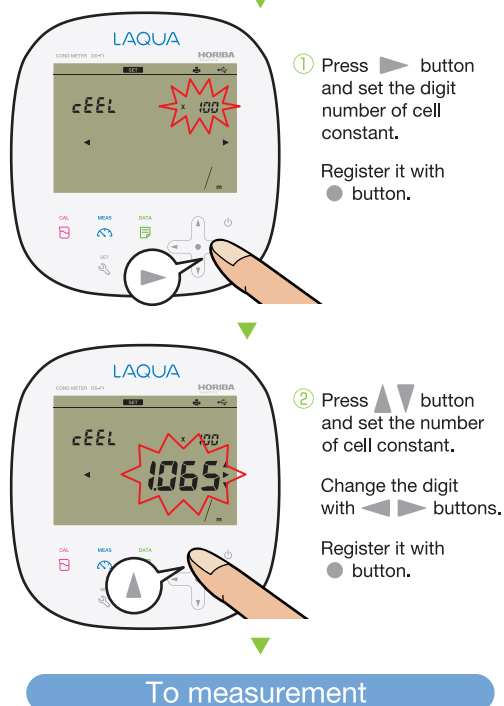

To measurement

Measurement

① Enter the measurement mode

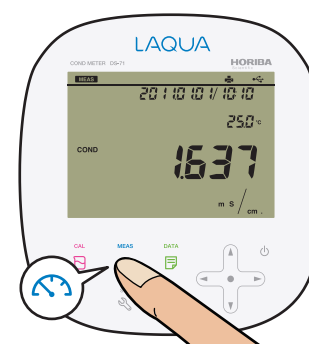

② Start the measurement



When the **HOLD** indication changes from the blinking state to the lit state, measurement is complete.